



1. Install small volume assembly into vessel plate
2. Insert correct small volume vessel into small volume assembly
3. Install small volume paddle or basket into dissolution machine
4. While the machine head is up drop one end of the DEPSET1015 tool, (10mm or 15mm) whichever your test calls for, into small volume vessel as shown in Figure 1
5. Lower the machine head so the paddle or basket touches the top of the DEPSET1015 tool
6. Lock your paddle or basket in place to ensure the height will remain the same
7. Raise the head and remove the DEPSET1015 tool
8. When ready to begin your test, your height will be set after lowering the head again